HP-B1 Basic Heating Plate

Please read the User Manual carefully before use, and follow all operating and safety instructions!

user manual

english

User Manual

HP-B1 Basic heating plate

Preface

Users should read this Manual carefully, follow the instructions and procedures, and beware of all the cautions when using this instrument.

Service

In order to guarantee this equipment works safely and efficiently, it must receive regular maintenance. In case of any faults, do not try to repair it yourself. If help is needed, you can always contact your supplier or Labbox via **www.labbox.com**.

Please provide the customer care representative with the following information:

- Serial number
- Description of problem
- Your contact information

Warranty

This instrument is warranted to be free from defects in materials and workmanship under normal use and service, for a period of 24 months from the date of invoice. The warranty is extended only to the original purchaser. It shall not apply to any product or parts which have been damaged on account of improper installation, improper connections, misuse, accident, or abnormal conditions of operation.

For claim under the warranty please contact your supplier.

Introduction

This product is suitable for use in various industries, agriculture, factories, medicine, and scientific research labs for heating. Other applications include household baking, boiling, and braising.

Characteristics

1. Designed in the latest style, it features stepless temperature control.
2. It is constructed with high-quality cold-rolled steel and has an electrostatically sprayed surface.
3. The fully enclosed heating plate offers the characteristics of flamelessness, rapid heating, excellent thermal insulation properties, a long service life, and reliability.

Main Technical Parameters

Voltage	220 V / 50 Hz
Heating Power	1500 W
Heating Plate Diameter	185 mm
Working Time	Continuous
Maximum temperature	$450 \pm 45^{\circ}\text{C}$

Operation Instructions

1. Connect the power cord to the machine, and the indicator light will illuminate.
2. Place the item that needs heating on the heating plate, then turn the temperature regulation knob clockwise to set the desired temperature. The heating plate will start working. A larger rotation range corresponds to a higher heating temperature.
3. Once the heating article reaches the required temperature, turn the knob in the opposite direction. Turn off the temperature regulation knob when it stops heating; this will stop the machine.

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